



## How to install the Sublicenses (from Web)

... Please turn page for information on how to deal with sublicenses and other files received on a USB Stick.

### Sublicenses for Analyzer 2


- 1) Register your dongle here: <http://www.brainproducts.com/productreg.php>
- 2) After having received your access data you can login and download your individual license-file for Analyzer 2 at <http://www.brainproducts.com/downloads.php> (at the very bottom of the page, see picture below)




- 3) Independently on your operating system (Windows XP, Windows Vista, Windows 7), install the license-file by double-clicking on the downloaded installation program (License\_zxxxxx.exe) and follow the wizard instruction. In the file name „z“ represents the type of dongle (U, UR, URA, N, VURA), „x“ represents digits.

### Sublicenses for Recorder (and Analyzer 1)

- 1) Register your dongle here: <http://www.brainproducts.com/productreg.php>
- 2) After having received your access data you can login and download the Recorder license-file at <http://www.brainproducts.com/downloads.php> (at the bottom of the page, see picture below)

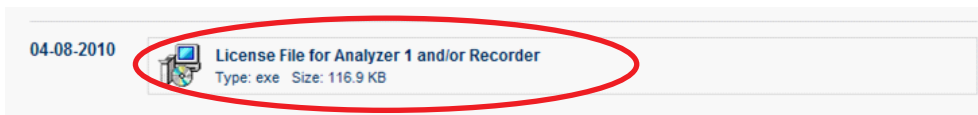

- 3) The installation depends on your operating system:
  - **Windows XP**  
When working with Windows XP, install the license-file by double-clicking on the downloaded installation program (BrainProducts-License-xxxxxxx-xxxx.exe) and follow the wizard instruction. In the file name „x“ represents digits.
  - **Windows Vista or Windows 7**  
When working with Windows Vista or Windows 7, click on the license-file icon with the right mouse button and select in the drop-down menu [Run as administrator] (see picture on the right).

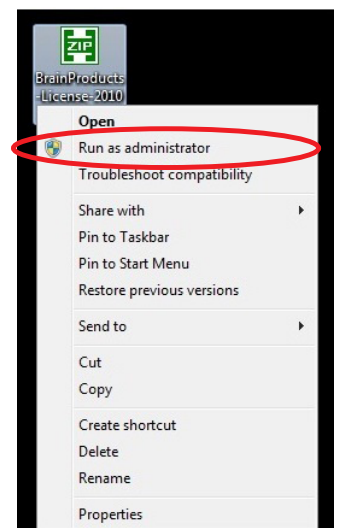



## How to install the Sublicenses (from USB Stick)

⚙️ Please turn page for information on how to deal with sublicenses downloaded from the internet

### Sublicenses for Analyzer 2

1) If you haven't done so yet, connect the dongle to your computer and install BrainVision Analyzer 2 and its updates from the enclosed BrainVision Installation DVD.

2) Connect the silver USB stick to your computer, unpack the file „License\_zxxxxx.zip“ and install the corresponding .exe file (License\_zxxxxx.exe). Make sure your dongle is connected when installing the file.

In the file name „z“ represents the type of your dongle (U, UR, URA, N, VURA), „x“ represents digits.

3) Independently on your operating system (Windows XP, Windows Vista, Windows 7), install the license-file by double-clicking on „License\_zxxxxx.exe“ and follow the wizard instruction.

### Sublicenses for Recorder (and Analyzer 1)


1) If you haven't done so yet, connect the dongle to your computer and install BrainVision Recorder and its updates from the enclosed BrainVision Installation DVD.

2) Connect the silver USB stick to your computer, unpack the file „BrainProducts-License-xxxxxx-xxxx.zip“ and install the corresponding .exe file (BrainProducts-License-xxxxxx-xxxx.exe).

In the file name „x“ represents digits.

3) The installation depends on your operating system:

- **Windows XP:** When working with Windows XP, install the license-file by double-clicking on „BrainProducts-License-xxxxxxx-xxxx.exe“ and follow the wizard instruction.  
In the file name „x“ represents digits.
- **Windows Vista or Windows 7:** When working with Windows Vista or Windows 7, click on the license-file icon with the right mouse button and select in the drop-down menu [Run as administrator] (see picture on the right).

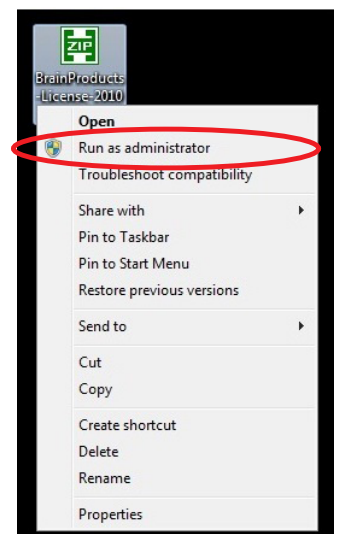

### Unlimited License

If your USB stick contains a file called „Update-zxxxxx-Unlimited.zip“, your dongle is currently time locked and will expire soon. Installing the file will remove the time lock.

1) Connect the silver USB stick to your computer, unpack the file „Update-zxxxxx-Unlimited.zip“ and install the corresponding .exe file (Update-zxxxxx-Unlimited.exe). Make sure your dongle is connected when installing the file.

In the file name „z“ represents the type of your dongle (U, UR, URA, N, VURA), „x“ represents digits.

2) Independently on your operating system (Windows XP, Windows Vista, Windows 7), install the file by double-clicking on „Update-zxxxxx-Unlimited.exe“ and follow the wizard instruction.

3) The installation depends on your operating system:

- **Windows XP:** When working with Windows XP, install the license-file by double-clicking on „BrainProducts-License-xxxxxxx-xxxx.exe“ and follow the wizard instruction. In the file name „x“ represents digits.
- **Windows Vista or Windows 7:** When working with Windows Vista or Windows 7, click on the license-file icon with the right mouse button and select in the drop-down menu [Run as administrator] (see picture above).