

## How to keep your PowerUnit fit and well

The PowerUnit is the new rechargeable lithium-ion battery from Brain Products. With an increased capacity and low self-discharge rate, it allows you significantly longer recording times and keeps its charge for a longer time in storage. The new PowerUnit is the standard battery for actiCHamp Plus, but it is also compatible with the actiCHamp.

Caring for your PowerUnit will ensure a reliable and long-lasting power source for your recordings. Be sure to follow some easy maintenance considerations to keep your PowerUnit fit and well:

- 1) Between scheduled recordings (within hours or days) the charger can stay connected to the battery.
- 2) If PowerUnit is not used for an extended period of time (weeks and months), disconnect the charger when the battery is fully charged.
- 3) To avoid deep discharge, fully recharge PowerUnit every four months.
- 4) During charging, operation and storage, keep the PowerUnit within a recommended temperature range of 10 °C to 30 °C (50 °F to 86 °F).
- 5) After use, disconnect PowerUnit from actiCHamp or actiCHamp Plus.
- 6) Only use the charger provided by Brain Products.