

Caution: To ensure correct and safe use of the products, make sure that you also read the Operating Instructions carefully prior to operation.



BRAINVISION
engineering

actiChamp-Plus Setup




- 1 Installing the software:** Insert the Brain Products Application Suite USB stick and install the **Recorder** software as described in the User Manual. The User Manual can be found on the Application Suite USB stick.

- 2 Installing the license dongle:** Plug in the USB dongle. The dongle driver is installed automatically. When installation is complete, the dongle will light up red.


Important: Your computer must be equipped with at least two USB2 ports for the software license dongle and the **actiChamp Plus** amplifier.




- 3 Optional step:** Insert the **actiChamp Plus** modules to upgrade your system. Use a suitable screw driver to undo and tighten the screws. Disconnect the **actiChamp Plus** completely before inserting modules.

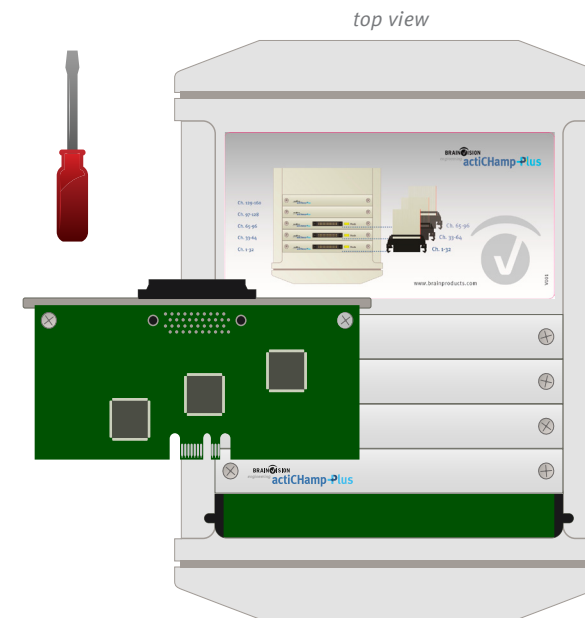

- 4** Connect the **actiChamp Plus** amplifier with the PowerUnit cable to the PowerUnit.

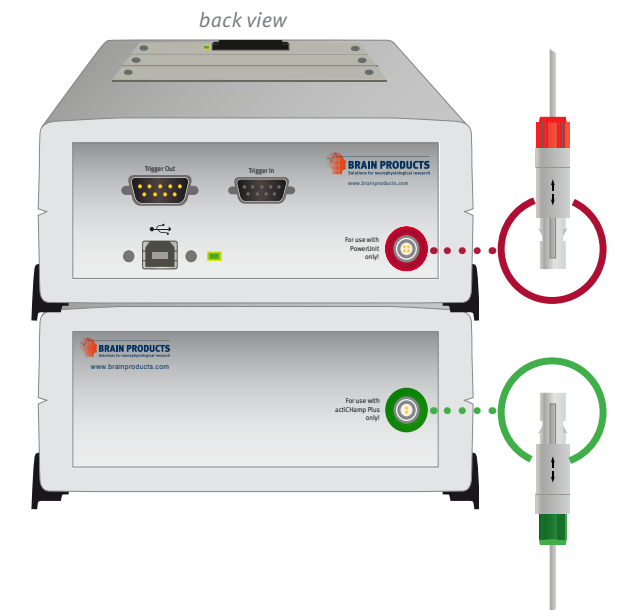


- 5** Connect the **actiChamp PI**

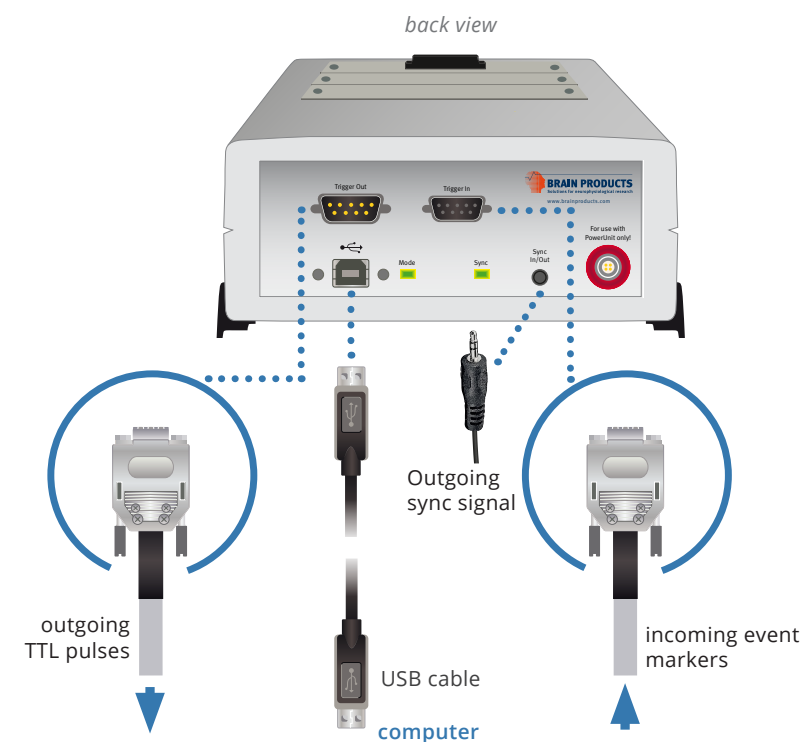


- 6** Connect the GND electrode to the **actiChamp Plus** amplifier.



- 7** Connect your electrodes to the **actiChamp Plus** amplifier.

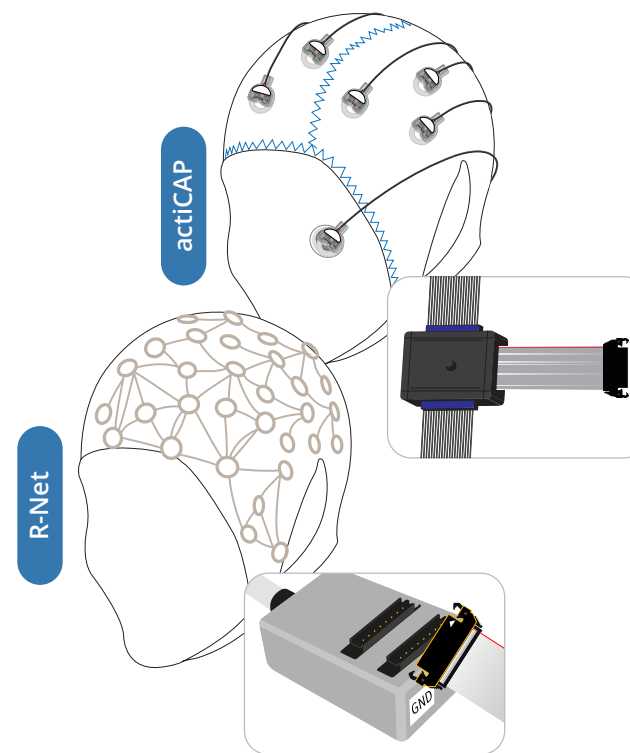


- 8** After use: Disconnect the PowerUnit from the **actiChamp Plus**. For information about correct battery handling, please refer to the **actiChamp Plus** operating instructions.

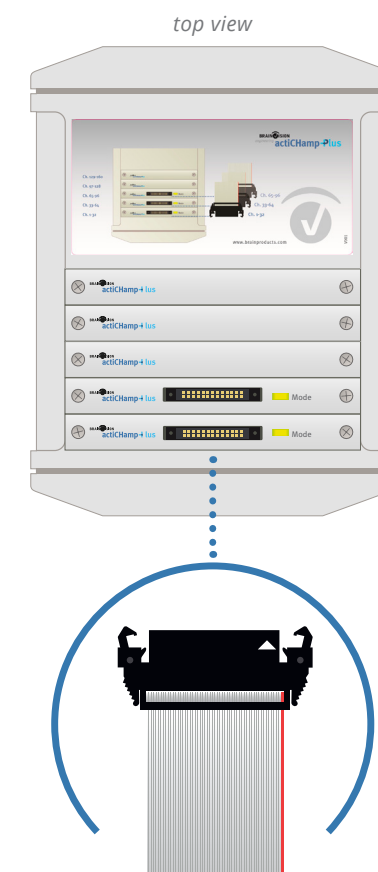

Setup Your actiChamp Plus System



Version 002
© Brain Products GmbH



Brain Products GmbH
Zeppelinstrasse 7
82205 Gilching
Germany
T +49 (0) 8105 733 84 0
F +49 (0) 8105 733 84 505
sales@brainproducts.com
www.brainproducts.com

Worldwide Caution:
Our products are scientific equipment for
INVESTIGATIONAL USE ONLY! Medical use
e.g. for diagnosis, treatment of disease or
other such purposes is strictly forbidden.