

Caution: To ensure correct and safe use of the products, make sure that you also read the Operating Instructions carefully prior to operation.



BRAINVISION professional **BRAINAMP Setup**




BRAIN PRODUCTS
Solutions for neurophysiological research

Brain Products GmbH
Phone +49 (0) 8105 733 84 0
sales@brainproducts.com
www.brainproducts.com

1

Installing the software: Insert the Brain Products Application Suite USB stick and install the **Recorder** software as described in the User Manual. The User Manual can be found on the Application Suite stick.

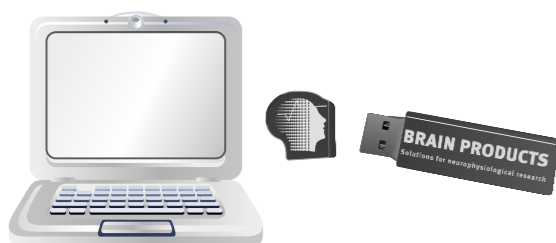

Plug in the USB stick and open the installation menu via double click on autorun.exe.

2

Installing the dongle: Plug in the USB dongle. The dongle driver is installed automatically. When installation is complete, the dongle will light up. Important: Your computer must be equipped with at least two USB2 ports for the software license dongle and the USB2 Adapter.





3

Installing the USB2 Adapter: Use the USB cable to connect the USB2 Adapter to your computer. Windows will detect the new hardware and install it automatically.




4

Installing the amplifier: Connect the amplifier to the USB2 Adapter using the fiber optic cable (FOC5 or FOC20). Important: Ensure that the connectors are plugged in correctly (amplifier 1 into fiber optic 1, amplifier 2 into fiber optic 2.)

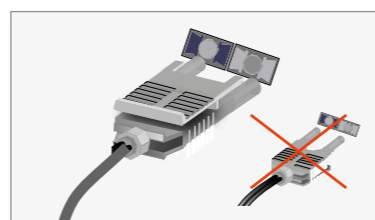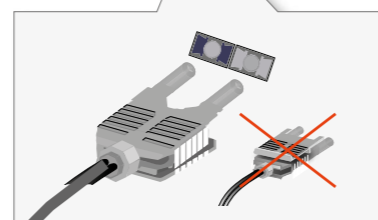

7

Connecting electrodes or electrode caps: Use the ribbon cable (BAC30 or BAC100) for connection.

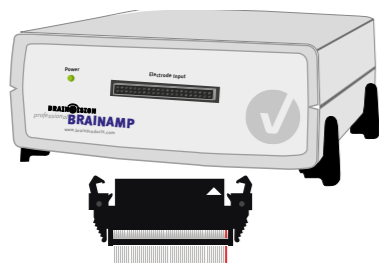


5

Connect the BrainAmp to the PowerPack using the PowerPack cable. Charge the PowerPack before use.



6

Switch on your amplifier (on the back). The power LED must light up.

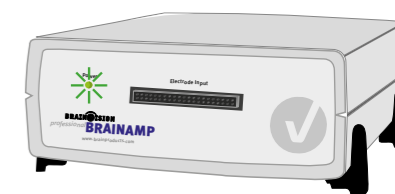


8

After use: Disconnect the PowerPack from the amplifier and immediately recharge the PowerPack. Leave the PowerPack connected to the charger if not in use. The PowerPack can be trickle charged and overcharging is not possible.

Setup Your BrainAmp System



© Brain Products GmbH



Brain Products GmbH
Zeppeinstrasse 7
82205 Gilching
Germany
T +49 (0) 8105 733 84 0
F +49 (0) 8105 733 84 505
sales@brainproducts.com
www.brainproducts.com

Worldwide Caution:
Our products are scientific equipment for
INVESTIGATIONAL USE ONLY! Medical use
e.g. for diagnosis, treatment of disease or
other such purposes is strictly forbidden.