



Setup Your LiveAmp Mobility Set





1 Check contents

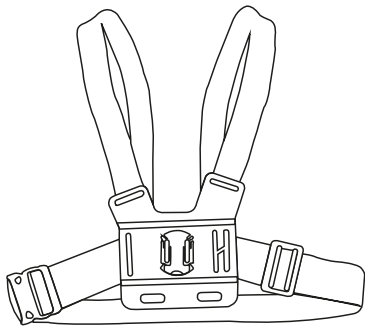


LiveAmp harness

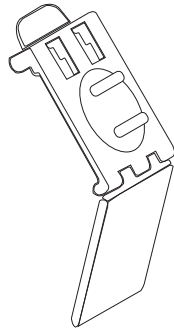

3x LiveAmp belt clips
Note: Belt clip may leave permanent marks on belt material

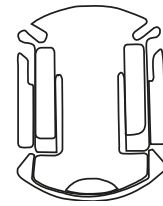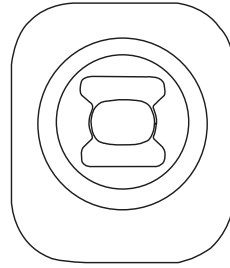


3x SPC mounts incl. microrail adapter

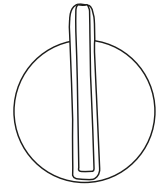


Stand tool to adjust the position



2 Prepare

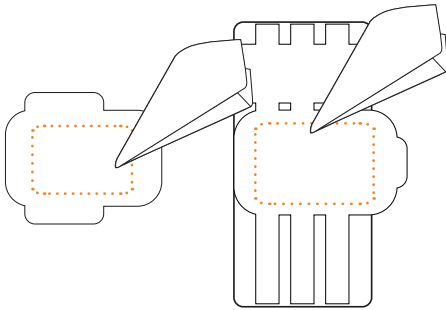

Clean the area where the SPC mounts go

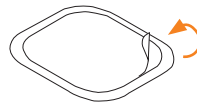


Stick the SPC mounts to the LiveAmp adapter

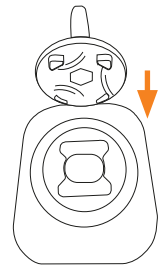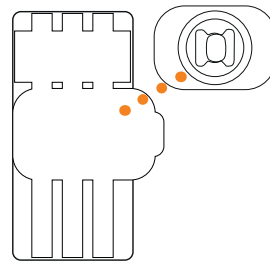


Note: Use stand tool to unlock the SPC mount, if necessary


3 Use

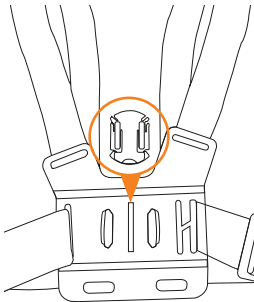


Place microrail adapter on LiveAmp harness

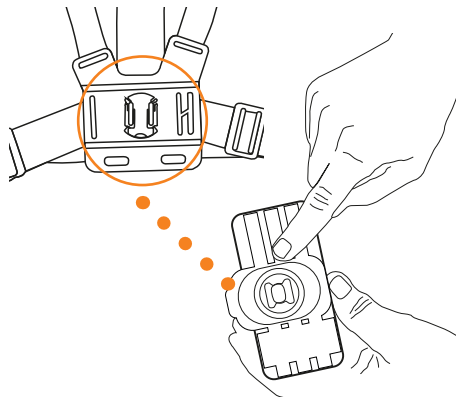

Place LiveAmp adapter with SPC mount onto the microrail adapter

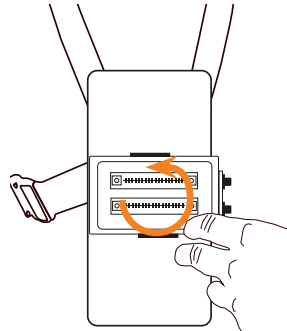

Turn 90° to secure the LiveAmp adapter

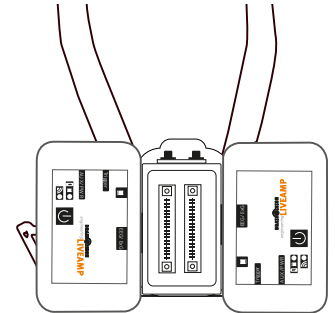

Exemplary LiveAmp 64 setup