

# TriggerBox Test I/O TriggerBox Driver



---

## Version 1.2.2

### **Windows® 10 driver update**

The driver has been updated for Windows® 10.

---

## Version 1.2.1

### **Event driven input port**

A new byte is available on the serial port only if the value of the TriggerBox input port has been changed. Instead of polling the input, the serial port event handling for received characters can now be used. The sampling interval of the input port is 5 ms, therefore to be recognized the trigger pulse width on the input port must be at least 5 ms.

---

## Version 1.1.1

### **Support for Windows® 10 and optimized support for MatLab®**

The new driver of the TriggerBox supports Windows® 10 and Matlab® as of 2015.  
This version is installed through the TriggerBox Test I/O software.

#### **Installing the new version**

1. Download the new **TriggerBox Test I/O 1.1.1** from:  
[www.brainproducts.com](http://www.brainproducts.com) > Downloads & Support > Downloads > Tools & Accessories.
  2. Provide your Brain Products user name and password.
  3. Click on the downloaded file to run the installation.
- ▶ Alternatively, you can run the installation of the **TriggerBox Test I/O 1.1.1** from the Application Suite DVD.

---

## Reporting errors and support

If you require support or if you discover a mistake in the manual, the software or during operation, please contact:

**Brain Products GmbH**  
**Zeppelinstraße 7**  
**D-82205 Gilching**  
**Phone: +49 8105 73384 – 0**  
**Fax: +49 8105 73384 – 505**  
**Web site: <http://www.brainproducts.com>**  
**Email: [techsup@brainproducts.com](mailto:techsup@brainproducts.com)**

On request, the support team will assist in clarifying faults and problems.

